

What's in the box?

Every matrIQ instrument will ship with the following in the box:

- The matrIQ instrument
- The matrIQ Quick Start User Guide
- A USB media drive containing the matrIQ installer and user documentation:
 - matrIQ Installer (**MatrIQ-1.X.X.exe**)
 - matrIQ Quick Start Guide
 - matrIQ instrument User Manual
- matrIQ AC/DC Power Adapter and cable
- USB-A to USB-B cable

The installer and user documentation can also be downloaded from:

www.coherent-solutions.com/resources/matrIQ/.

Getting Started




IMPORTANT

If **another matrIQ instrument is already connected** to the client computer **over USB**, consult the User Manual on how to **change the USB IP address**.


matrIQ can be communicated with over **Ethernet** or **USB** connection. Both connection methods will allow control of the instrument through the cohesionUI™ graphical user interface and with SCPI commands.

STEP ONE – Power on the instrument

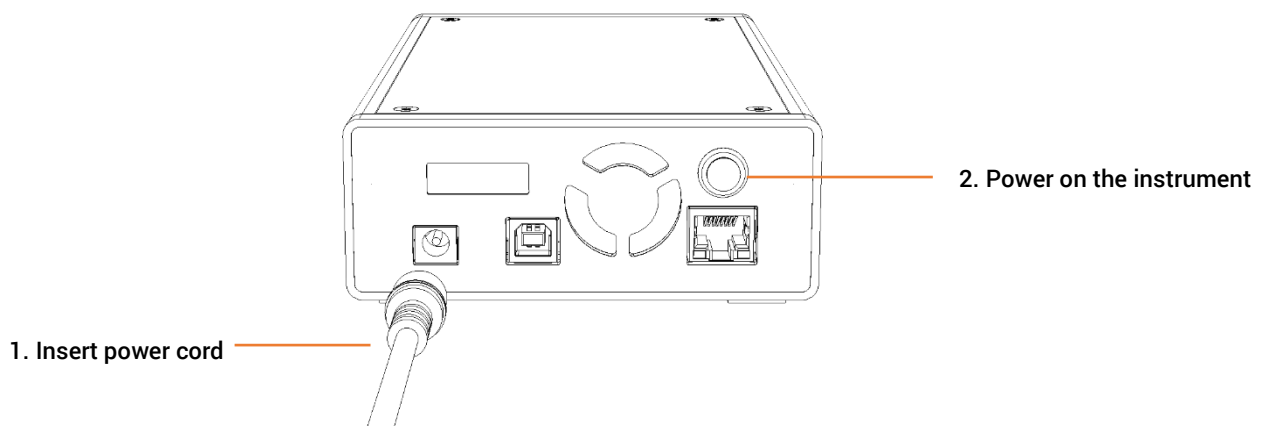


STEP TWO

Option 1: Communication via USB

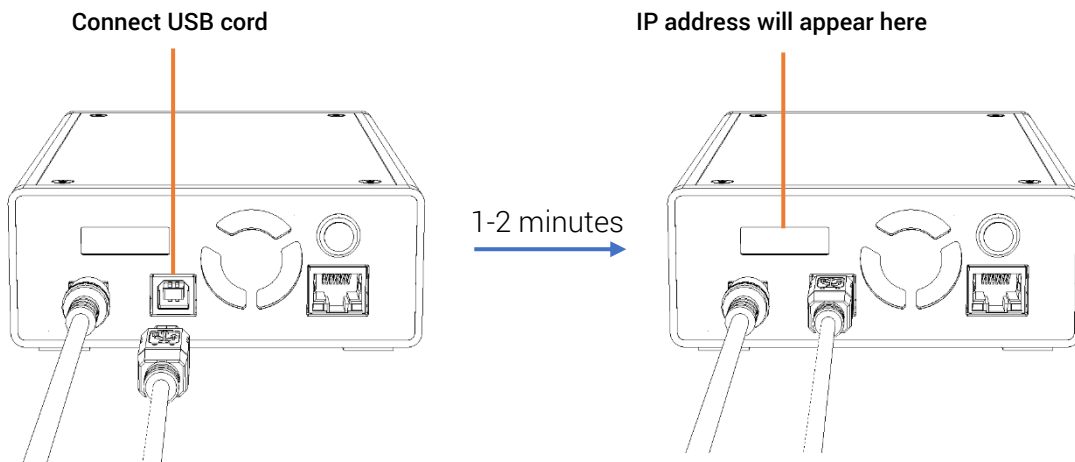

Option 2: Communication via Ethernet

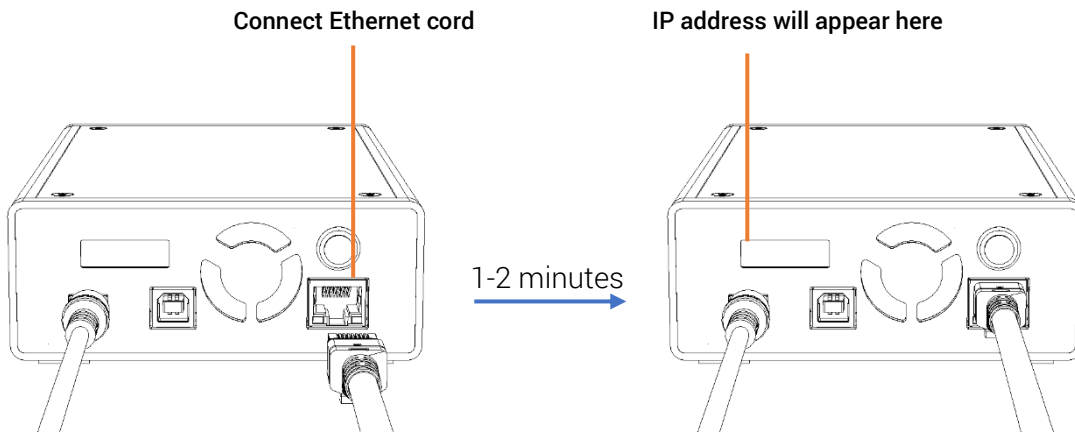

STEP THREE – Run the matrIQ Installer

Double click and run the **MatrIQ-1.X.X.exe** matrIQ Installer from the provided USB Media drive.

STEP FOUR – Run the matrIQ application

Double click and run the **MatrIQ** desktop application.



Follow the on-screen prompts from the MatrIQ application landing page to use cohesionUI to control the matrIQ instrument.



NOTE

Consult the provided matrIQ instrument User Manual for further information on configuring the Ethernet/USB IP address.